

The background is a dark purple gradient. On the left, there are two vertical neon lines, one blue and one magenta, with a small horizontal magenta line intersecting the blue one. On the right, a large, glowing magenta arc curves from the top towards the bottom. The text 'AVEVA WORLD' is centered in a white, bold, sans-serif font.

AVEVA WORLD

APRIL 9, 2025

---

## Cytiva & SyVento: Enhancing Biomanufacturing Efficiency with FlexFactory™ Historian

SESS-146

Tim Wortley – Cytiva



# Enhancing Biomanufacturing Efficiency with FlexFactory™ Historian

**A Case Study from SyVento Biotech**

Tim Wortley, Cytiva

# Who are we



## Vision

Our vision is a world in which access to life-changing therapies transforms human health.

## Mission

Our mission is to advance and accelerate therapeutics.



## Encapsulating your ideas

SyVento BioTech is a biotechnology company offering comprehensive services in the development of lipid-based nanocarriers and LNP technology. We specialize in the development of lipid-based formulations for pharmaceutical products, including RNA-based drug products.

# Why it matters ... Let's meet Emily







**FOCUS  
FORWARD**

**SHORT FILMS / BIG IDEAS**

# Introduction

Cytiva's FlexFactory™ historian biomanufacturing solution built on top of AVEVA™ PI System transforms mRNA manufacturing at Syvento™ Biotech in Poland.

A discussion on the challenges in mRNA manufacturing and the transformative improvement on efficiency, costs and regulatory compliance from data management



# The regulatory landscape and market trends

## Why comply?

- Reputation, costs (fines), stay in business, access to global markets, directors & employees stay out of jail...
- Well recognized guidance documents on how to deliver safe and effective drugs. Similar standards apply globally.

## How to comply (data)

- “If it didn’t get documented (reviewed, approved, signed), it didn’t happen”
- The (process, instrument, analytical, manufacturing) **data** is sacrosanct, must be original and proven tamperproof.

## Complex Trends

- “Zoo” of biotherapeutics (mAbs, mRNA, AAV, pDNA, siRNA, EBT)
- Faster, cheaper, simpler (Viz SARS-COV2 vaccine in ~11months<sup>2</sup>, typically many years<sup>1,3</sup>)



1. [The development of COVID-19 vaccines in the United States: Why and how so fast? – PMC](#)
2. Emergency authorization Dec 2020, Approval Dec 2021 [COMIRNATY | FDA](#)
3. Merck’s Gardasil was in development at U-Queensland in 1991 to approval in 2014 (23 years), [Gardasil -Wikipedia](#)



# Bioprocessing is challenged with inefficiency

Over 300 error-prone manual interventions in average 2000L mAb SU process

More than 80% of company time is spent collecting and cleaning data sets<sup>1</sup>

10–20% of the time biopharma data can't be accessed<sup>2</sup>

Success of initial production run can be as low as 47%<sup>3</sup>

Biopharma scores lower than other industries in adoption of automation<sup>4</sup>

High batch failure rates and manufacturing deviations due to human error<sup>5</sup>



<sup>1</sup> <https://www.outsourcing-pharma.com/Article/2019/01/09/AI-and-improvements-to-pharma-manufacturing#>

<sup>2</sup> [https://www.idbs.com/wp-content/uploads/2020/12/IDBS\\_Agility\\_Article\\_Sept\\_2020\\_V4.pdf](https://www.idbs.com/wp-content/uploads/2020/12/IDBS_Agility_Article_Sept_2020_V4.pdf)

<sup>3</sup> <https://www.biopharminternational.com/view/review-exception-connecting-dots-faster-batch-release>

<sup>4</sup> <https://www.mckinsey.com/industries/pharmaceuticals-and-medical-products/our-insights/automation-and-the-future-of-work-in-the-us-biopharma-industry>

<sup>5</sup> Bioplan Associates, "17th Annual Report and Survey of Biopharmaceutical Manufacturing Capacity and Production," April 2020: [https://bioplanassociates.com/wp-content/uploads/2020/06/17th-Annual-Biomfg-Report\\_TABLE-OF-CONTENTS-2020.pdf](https://bioplanassociates.com/wp-content/uploads/2020/06/17th-Annual-Biomfg-Report_TABLE-OF-CONTENTS-2020.pdf)

# Automation and control can lead to large efficiency gains

**Up to 75%**


reduction in paper handling and storage<sup>1</sup>

**Up to 20%**

increase in employee efficiency<sup>3</sup>

**Up to 38%**

reduction in labor<sup>5</sup>

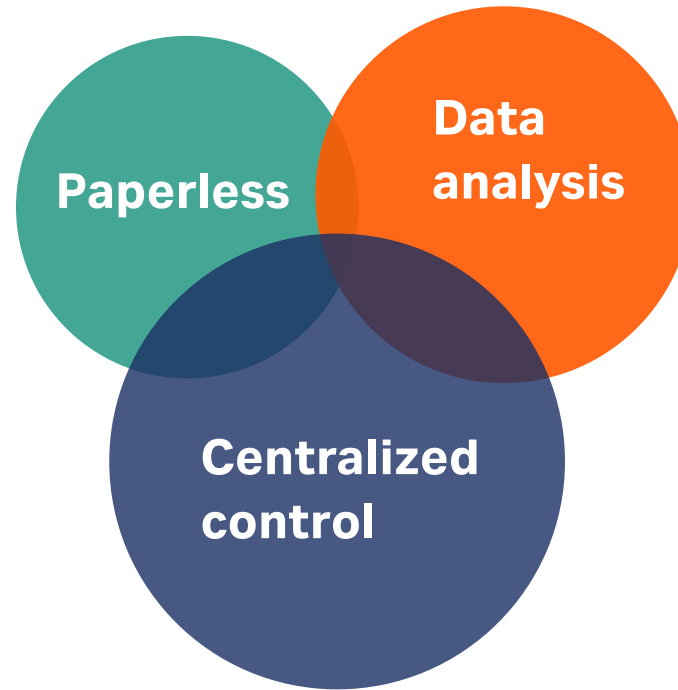

**Up to 85%**

reduction in batch review time<sup>2</sup>

**Up to 79%**

reduction in failed batches<sup>4</sup>

**Up to 67%**

reduction in deviations<sup>5</sup>

**Enable electronic data capture, storage, access, and analysis**

1, 2. [http://www.sysiss.com/fileupload/sysiss\\_documents/Systematic-Takeda%20EBRS%20case%20studyV1-600dpi.pdf](http://www.sysiss.com/fileupload/sysiss_documents/Systematic-Takeda%20EBRS%20case%20studyV1-600dpi.pdf)

3. <https://www.mckinsey.com/industries/pharmaceuticals-and-medical-products/our-insights/automation-and-the-future-of-work-in-the-us-biopharma-industry#>

4. Bioplan Associates, "17th Annual Report and Survey of Biopharmaceutical Manufacturing Capacity and Production," April 2020: [https://bioplanassociates.com/wp-content/uploads/2020/06/17th-Annual-Biomfg-Report\\_TABLE-OF-CONTENTS-2020.pdf](https://bioplanassociates.com/wp-content/uploads/2020/06/17th-Annual-Biomfg-Report_TABLE-OF-CONTENTS-2020.pdf)

5. Cytiva – Biopharm Services study using BioSolve™, unpublished, March 2022

# mRNA therapeutics

Synthetic messenger RNA molecules are introduced and temporarily (hours-days) hijack our cells function to build the programmed target<sup>1</sup> to extend our immune system e.g.

- Infectious diseases (e.g. MERS-CoV, SARS-CoV2)
- Anti-cancer molecules / Oncology vaccines
- Cell & Gene therapies (CRISPR, CAR-T)
- Other vaccines (Rabies, HIV, Nipah etc)

## Advantages

- **Rapid Development:** mRNA-based treatments can be designed and produced quickly, enabling swift responses to emerging health threats, such as pandemics<sup>2</sup>
- **Efficient Production:** mRNA therapeutics are produced in cell-free systems, which simplifies manufacturing and reduces resource intensity compared to traditional protein-based biologics<sup>3</sup>
- **Versatility:** mRNA can encode a wide range of proteins, making it adaptable for various applications, including vaccines, cancer immunotherapy, and protein replacement therapies<sup>4</sup>

## Pandemic response

10–100 L IVT scale  
<1–300 kg/year



## Other infectious diseases and oncology

0.1–10 L IVT scale  
<1 g–50 kg/year



## Personalized therapies

10–50 mL IVT scale  
<1 g/year



# Future proofing global manufacturing

## flexible solutions required in-region-for-region

- mRNA technologies are relatively new (BioNTech founded 2008)
- Obscure (lost in Zoo?) until Covid
- Pandemic demonstrated need for in-region-for-region vaccine manufacturing
- mRNA vaccines can be mass-produced in a trailer.

### Considerations

- Different process needs with different mRNA types
- No (new) platform process
- Rapidly evolving technology
- Unique problems need novel solutions (e.g. lipid nanoparticles)
- Few approved therapies
- Capital and operating costs critical.
- Substantial differences between pandemic response versus personalized therapeutics




# mRNA process – unique challenges

## Process intensification

- 1-2 day process, many small batches (typical mAB is 15-30day cycle)
- Documentation, batch release (300 batches/year versus 30)
- High yield processes – 1L reactor  $\rightarrow$  4g mRNA<sup>2</sup>  $\rightarrow$  100,000 x 30 $\mu$ g dose<sup>1</sup>

## Process needs

- Process, consumables, reagents, data usage 10-100x faster than conventional
- Paper processes (and staff) can no longer keep up, especially in a high stress (pandemic?) environment
- 2 day manufacturing process waiting for 2 month batch release not acceptable when patients are dying.
- Release by Exception

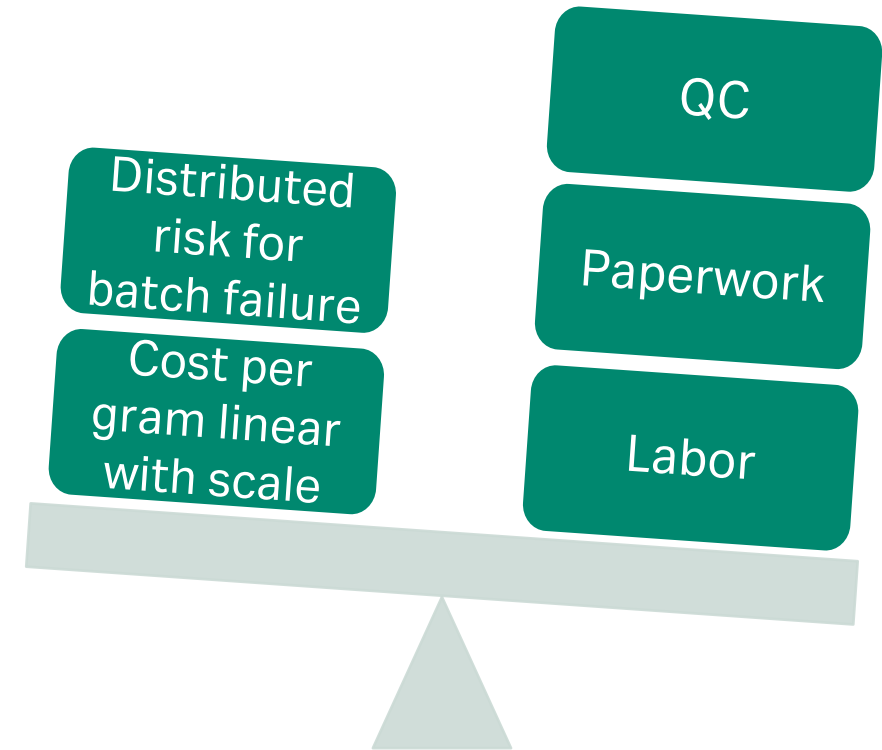




# Intensified and complex manufacturing

The mRNA future manufacturing will have to handle multiple products and multivalent products.


Challenge:

- short process and intensified manufacturing
- line clearance and batch-to-batch separation
- maximize facility utilization

mRNA multiproduct scheduling example

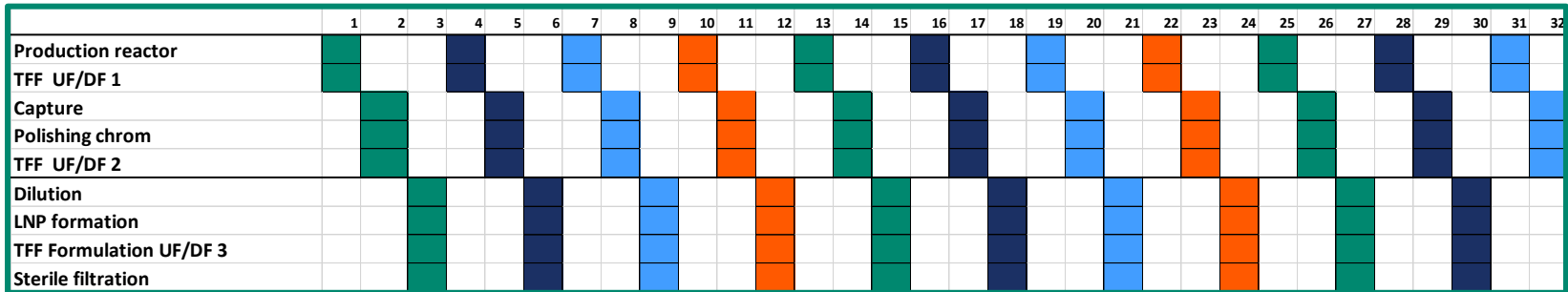

mRNA multivalent product scheduling example

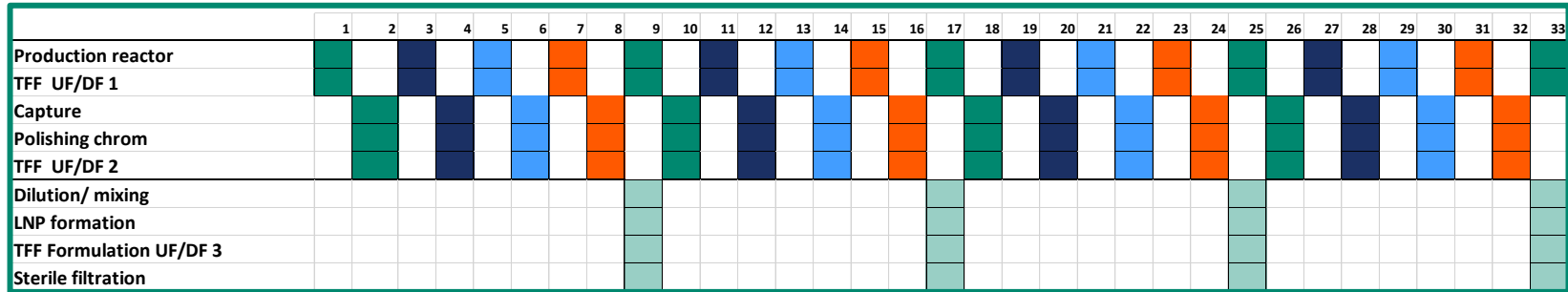

# How historian technologies and pre-configured, pre-verified solutions improve manufacturing outcomes

## Standard Cytiva and third party equipment

- Custom designed equipment is typical in industry, and not necessary
- Modern equipment have standard interfaces that are accommodated by PI Connectors and Interfaces2

## Pre-design, pre-verified and documented solution

- Client receives an integrated solution from a single vendor that can rapidly configure, deploy and verify an industry standard expandable and flexible solution


## Standard, templated and automated AVEVA PI System build.

- PI System support can be deployed for standard equipment in hours, obviating need for typical design-specification-build-test-document life cycle
- Standard PI System allows customers and their integrators to modify, extend and expand the deployed system.

## Comprehensive CQV<sup>1</sup> documentation, protocols and support

- Working with standard equipment and software configuration optimizes the CQV process.
- Typically double+ the automation cost -- reduces significantly

Cytiva

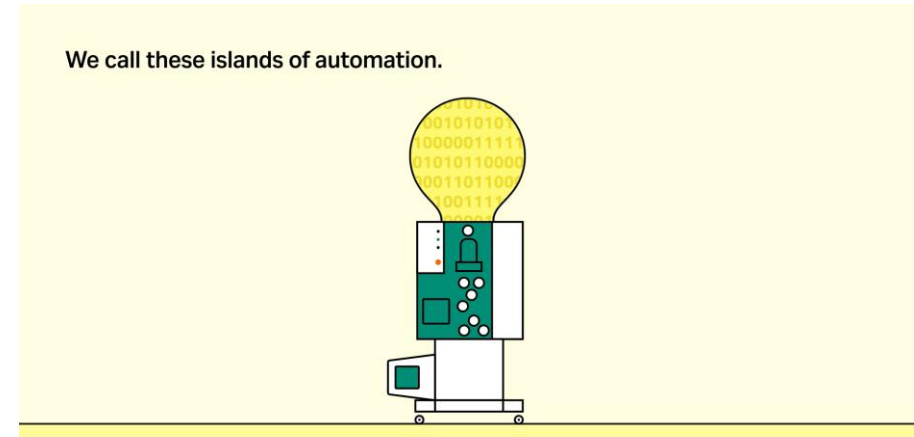


1. [Explaining CQV, C&Q, and Commissioning and Qualification | Kneat](#)

# Typical FlexFactory historian deliverables




## Specifications

- General (Hardware)
- Functional

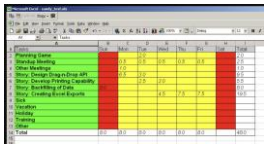

## Configuration Specification (Template)

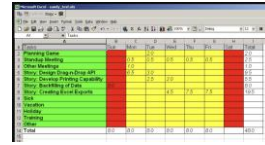

## Configuration Specification

- TOP documentation
- Semi-Auto generated
- Tag & AF lists
- Test results docs
- Configuration docs

Automatically

generated




## Standard Build/ Work instructions


Process Equipment Config  
PI System configuration

- Tags
- Audit trail interface
- Asset Framework

Cytiva

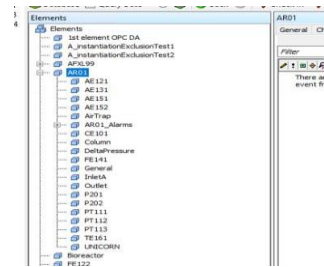


## Tags, connectors and Interfaces

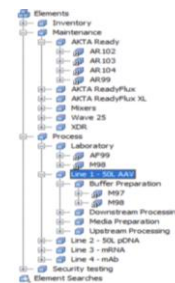


## PIAF

## Infrastructure Build

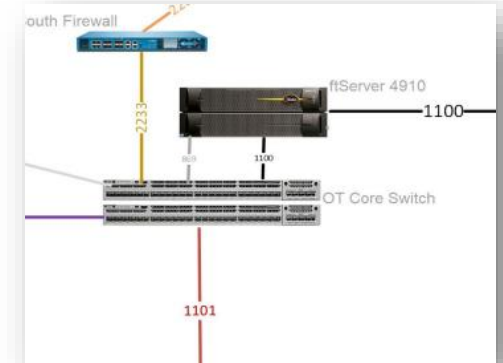

## PI Vision Displays

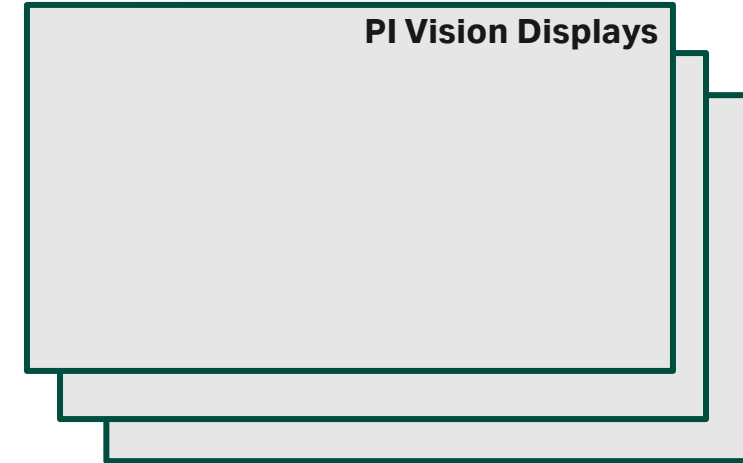

## Turn-over-Package (Commissioning & Qualification documentation)



# Conclusion

## **FlexFactory historian powered by AVEVA PI System has delivered world class mRNA manufacturing capacity to Poland**

- Syvento can deliver life saving therapeutics of a wide range of types, ensuring advanced treatments reach patients who need them.
- Improve access to personalized medicines and mRNA therapies across Europe to meet future CAR-T, oncology and pandemic vaccine needs.
- Enable creation of a new mRNA center of excellence, supporting global therapeutic drug research, development and manufacturing.

## **The FlexFactory provides a future-proofed and digitally enabled manufacturing platform**

- Meet future regulators expectations for data integrity and analysis.
- Provide frictionless movement of data between manufactures, client biopharmaceuticals and regulators
- Provide for more rapid and smarter drug batch release, improve therapeutic drug development time to market.





Cytiva and the Drop logo are trademarks of Global Life Sciences IP Holdco LLC or an affiliate.  
FlexFactory and AKTA, are trademarks of Global Life Sciences Solutions USA LLC or an affiliate doing business as Cytiva.

Syvento is a trademark of Spółka SyVento Sp. All other third-party trademarks are the property of their respective owners.

© 2025 Cytiva

All goods and services are sold subject to the terms and conditions of sale of the supplying company operating within the Cytiva business. A copy of those terms and conditions is available on request.  
Contact your local Cytiva representative for the most current information.

For local office contact information, visit [cytiva.com/contact](https://www.cytiva.com/contact)



# Thank you

**Firstname Lastname**

emailaddress@cytiva.com

0.000.000.0000